

Stock Solution



TD-S Revision 2.0

Creation Date: 4/8/2019
Revision Date: 10/21/2019

15 mg/ml Luciferin Stock Solution for *In Vitro* Use (D-Luciferin & L-Luciferin)

Instructions

1. For a 15 mg/ml solution, dissolve 1.0 grams of D-Luciferin* into 66.6 ml sterile water.
2. Aliquot into tubes.
3. Freeze solution at -80°C (Good for up to 1 year).

Note: *Applies to D-Luciferin Potassium Salt, D-Luciferin Sodium Salt and L-Luciferin Potassium Salt (GoldBio Catalog # [LUCK](#)[®], # [LUCNA](#) or # [L-127](#)) for *in vitro* use.