

Stock Solution



TD-S Revision 2.1

Creation Date: 6/26/2014

Revision Date: 7/14/2022

100 mg/ml Ampicillin Stock Solution

Instructions

1. Weigh 1 g of Ampicillin ([Ampicillin \(Sodium\), GoldBio Catalog # A-301](#) [CAS 69-52-3, mw. = 371.4]).
2. Add 10 ml of sterile H₂O. Dissolve completely.
3. Prewet a 0.22 µm syringe filter by drawing through 5-10 ml of sterile H₂O and discard water.
4. Sterilize Ampicillin Stock through the prepared 0.22 µm syringe filter.
5. Stock may be kept at -80°C for 1 year.

Note: Ampicillin solutions degrade 13% after one week at -20°C.

Gold Biotechnology

St. Louis, MO

Ph: (800) 248-7609

Web: www.goldbio.com

Email: contactgoldbio86@goldbio.com