

Stock Solution



TD-S Revision 2.0

Creation Date: 4/8/2019
Revision Date: 10/17/2019

Coelenterazine Stock Solution for *In Vivo* Use

Instructions

1. Add 5 mg coelenterazine ([GoldBio Catalog # CZ](#)) to 1 ml of acidified methanol (add 50 μ l of concentrated HCl to 10 ml methanol).
2. Aliquot and store at -80°C for up to 4 weeks.
3. For use, thaw tube and add 20 μ l of coelenterazine stock solution to 130 μ l DPBS and inject immediately.