


## CY5-NHS Ester Technical Information

Product No.: B-430

Product Name: [CY5-NHS Ester](#)

Chemical Structure:

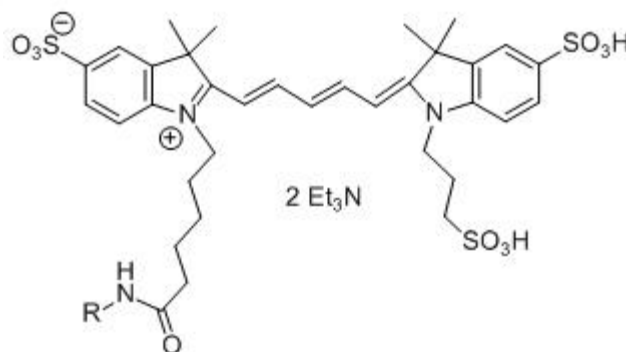

Chemical Composition: C<sub>38</sub>H<sub>46</sub>N<sub>3</sub>O<sub>11</sub>S<sub>3</sub>

Molecular Weight: 1050.97 g/mol (triethylammonium salt)

Solubility: Water, DMF, DMSO

Spectral Properties:

Absorbance/Emissions	649/670 nm
$\epsilon_{\max}$	250,000 M <sup>-1</sup> cm <sup>-1</sup>
CF <sub>280</sub>	0.05
CF <sub>260</sub>	0.05

Storage: Store at -20°C. (Product shipped at ambient temperature.)

### Introduction

Cy5-NHS Ester is a bright, photostable, and pH insensitive fluorescent dye that enables simple and efficient Cy5 fluorescent labeling of any primary amine-containing macromolecule. They are excited to about 98% of maximum with a krypton/argon laser (647 nm line) or to about 63% of maximum with a helium/neon laser (633 nm line). NHS activated esters react efficiently with primary amino groups (-NH<sub>2</sub>) forming stable amide bonds. When labeling antibodies or other proteins with Cy dyes, the brightest conjugates have dye/protein ratios ranging from 4-12<sup>(1)</sup>, depending on the specific applications. High ratios can increase non-specific background, loss of antibody binding affinity, and/or lead to aggregation. For this reason, optimal labeling ratios should be experimentally verified for any given application.

### Gold Biotechnology

St. Louis, MO

Web: [www.goldbio.com](http://www.goldbio.com)

Ph: (314) 890-8778

email: [contactgoldbio86@goldbio.com](mailto:contactgoldbio86@goldbio.com)

### Important Product Information

- NHS esters are moisture-sensitive. To avoid moisture condensation onto the product always let vial come to room temperature before opening; be careful to limit exposure to. The NHS-ester moiety readily hydrolyzes and becomes non-reactive; therefore, prepare stock solutions immediately before use. Prepare stock solutions in anhydrous solvents (e.g. DMSO or DMF) immediately before use.
- Hydrolysis of the NHS ester is a competing reaction in aqueous buffers. Conjugation with primary amines of proteins/peptides (i.e., acylation) is favored in the pH range 7-9 with concentrated protein solutions (e.g. 1-10 mg/ml). Use non-amine-containing buffers at pH 7-9, e.g. 100mM sodium phosphate, 150mM sodium chloride, pH 7.5, 100mM HEPES, pH 7.5, 100mM carbonate/bicarbonate, pH 8-9 or 100mM borate buffer, pH 8-9.
- Do not use buffers that contain primary amines, (e.g., Tris, glycine).

### Additional Materials Required

- Water-miscible organic solvent such as dimethyl sulfoxide (DMSO) or dimethyl formamide (DMF)
- Reaction buffer: Sodium phosphate buffer (100mM sodium phosphate, 150mM sodium chloride, pH 7.5).
- Desalting Spin Columns or Gravity Flow Desalting SEC Columns (e.g. Sephadex G-25).

### Procedure for Labeling an Antibody with Cy5-NH<sub>2</sub>

1. Dissolve antibody into a suitable amine-free buffer (e.g. 100mM sodium phosphate, 150mM sodium chloride, pH 7.5) to obtain a 1-5 mg/ml solution. If necessary buffer exchange/desalt into 100mM sodium phosphate, 150mM sodium chloride, pH 7.5 prior to labeling if residual traces of Tris or glycine contaminants are present.
2. Immediately before use, dissolve Cy5-NHS Ester reagent into a suitable volume of DMSO or aqueous buffer in order to achieve a reagent concentration 400-fold greater than the initial concentration of antibody. For example, if the IgG to be labeled is at 1 mg/ml (i.e. 6.7 $\mu$ M) dissolve the Cy5-NHS Ester reagent at 2.7mM in DMSO, if IgG is at 2.5 mg/ml dissolve the reagent at 6.75mM, etc.
3. Add the appropriate volume of NHS reagent to the antibody in order to achieve the desired molar excess of Cy5-NHS Ester reagent. To target a specific degree of labeling, refer to Figure 1 under calculations.
4. Incubate the reaction at room temperature for 60 minutes (or longer).
5. Remove excess non-reactive reagent by dialysis or desalting spin column into buffer of choice.
6. Measure absorbance of Cy5-antibody conjugate at 280 nm and 550 nm.
7. Determine Cy5 DOL and conjugate concentration. See Calculations.

### Gold Biotechnology

St. Louis, MO

Web: [www.goldbio.com](http://www.goldbio.com)

Ph: (314) 890-8778

email: [contactgoldbio86@goldbio.com](mailto:contactgoldbio86@goldbio.com)

### Calculations

Calculate degree of labeling (DOL) and Cy5 conjugate concentration (mg/ml) with the following Equations:

$$\text{Eq. 1} \quad \text{number of Cy5 dye per protein} = \frac{\text{molarity Cy5 dye}}{\text{molarity protein}}$$

$$\text{Eq. 2} \quad [\text{molarity of Cy5 dye}] = \frac{A_{650}}{\epsilon_{650}}$$

$$\text{Eq. 3} \quad [\text{molarity of protein}] = \frac{A_{280C}}{\epsilon_{280}}$$

$$\text{Eq. 4} \quad \text{mg/mL} = \frac{A_{280} - (650 \times 0.05)}{\left(\frac{E1\%}{10}\right)} \times \text{dilution factor}$$

$A_{550}$  = conjugate absorbance at 650 ± 2 nm

$\epsilon_{550}$  = molar extinction coefficient Cy5 dye = 250,000 M<sup>-1</sup> cm<sup>-1</sup>

$A_{280}$  = conjugate absorbance at 280 nm

$A_{280C}$  = corrected conjugate absorbance at 280 nm =  $A_{280} - (A_{650} \times (0.05))$

$\epsilon_{280}$  = molar extinction coefficient protein (M<sup>-1</sup>cm<sup>-1</sup>) =  $\frac{MW_p \times E1\%}{10}$

Example 1: Determine Cy5 DOL and conjugate concentrations as follows:

A Goat IgG antibody 0.1 mL at 1.0 mg/mL was labeled using a 20-fold molar excess Cy5-NHS reagent. After removal of excess dye with a spin column, the conjugate's A280 and A650 were determined in PBS (1:5 dilution), A280 = 0.2680 and A650 = 1.3783. Goat IgG M.W. 150kDa, E1% = 13.6 or 204,000 M<sup>-1</sup>cm<sup>-1</sup>).

**Calculate DOL (# Cy 5/protein) as follows:**


**By Equation 2**      molarity of Cy5 =  $\frac{1.3783}{250,000 \text{ M}^{-1}\text{cm}^{-1}}$  = 5.51 μM

**By Equation 3**      molarity of IgG =  $\frac{0.2680 - (1.3783 \times 0.05)}{204,000 \text{ M}^{-1} \text{ cm}^{-1}}$  = 0.877 μM

**By Equation 1**      Number of Cy5 per IgG =  $\frac{5.51 \mu\text{M}}{0.877 \mu\text{M}}$  = 6.28

**Calculate Cy5 -IgG protein concentration (mg/mL)**

**By Equation 4**      mg/mL =  $\frac{0.2680 - (1.3783 \times 0.05)}{1.36} \times 5$  = 0.066 mg/mL

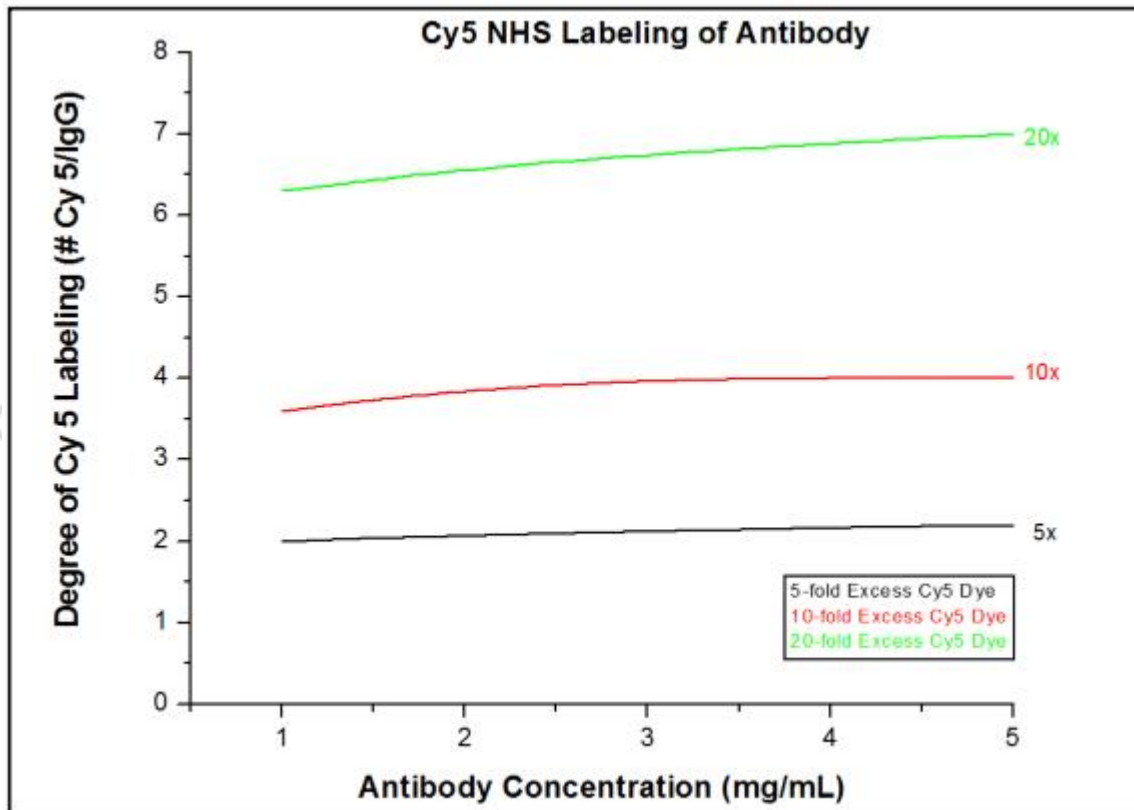

**Figure 1.** Goat IgG samples (100mM sodium phosphate, 150mM sodium chloride, pH 7.5) at various concentrations were labeled with 5, 10, or 20-fold excess Cy5-NHS Ester for 1 hour at room temp. Resulting conjugates generated the curves seen above. These can be used to target other dye levels under the buffer, pH, and solvent (DMSO) conditions used in this protocol.

## References

1. Southwick, P.L., et al. (1990). Cyanine dye labeling reagents-carboxymethylindocyanine succinimidyl esters. *Cytometry*. 11(3): 418-430.

*Cy is a trademark of GE Healthcare.*

**For research use only**

**Gold Biotechnology**

St. Louis, MO

Web: [www.goldbio.com](http://www.goldbio.com)

Ph: (314) 890-8778

email: [contactgoldbio86@goldbio.com](mailto:contactgoldbio86@goldbio.com)